

特色

美國原裝進口

採用 Quick-Connect 設計，安裝超快速

含調壓閥

Benefits

**Benefits** of using the **Autofil™** filtration system

- Vacuum hose is connected just one time
- Quick-Connect™ pedestal offers fast connection and hands-free operation
- Forward facing graduation marks on tubes and bottles
- Vacuum on and off switch conveniently located
- **Autofil™** design including an ergonomic bottle shape, molded finger grips and patented bottle cap
- Conveniently sold presterilized
- **Autofil™** is available as a complete system or as a bottle-top filter

**A** • Centrifuge tubes available in 15ml and 50ml  
• 0.1 µm, 0.22 µm and 0.45 µm PES membrane pore-size options available

**B** • Bottles available in 250ml, 500ml and 1000ml sizes  
• Available in 0.1 µm, 0.22 µm or 0.45 µm asymmetric membrane filter



**Easy initial setup of Autofil™ filtration system**



**1** Connect a vacuum tube just once to the back of the Quick-Connect™ pedestal.

**2** Vacuum connects automatically when filter is placed on the Quick-Connect™ pedestal.

**3** Once fluid is added to the top of the **Autofil™** system, turn the vacuum switch on.